

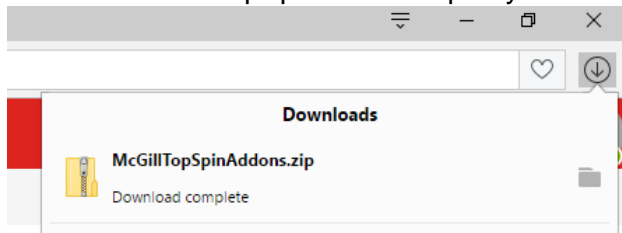



Installing McGill TopSpin Addons

1. Save the file McGillTopSpinAddons.zip to your Downloads directory

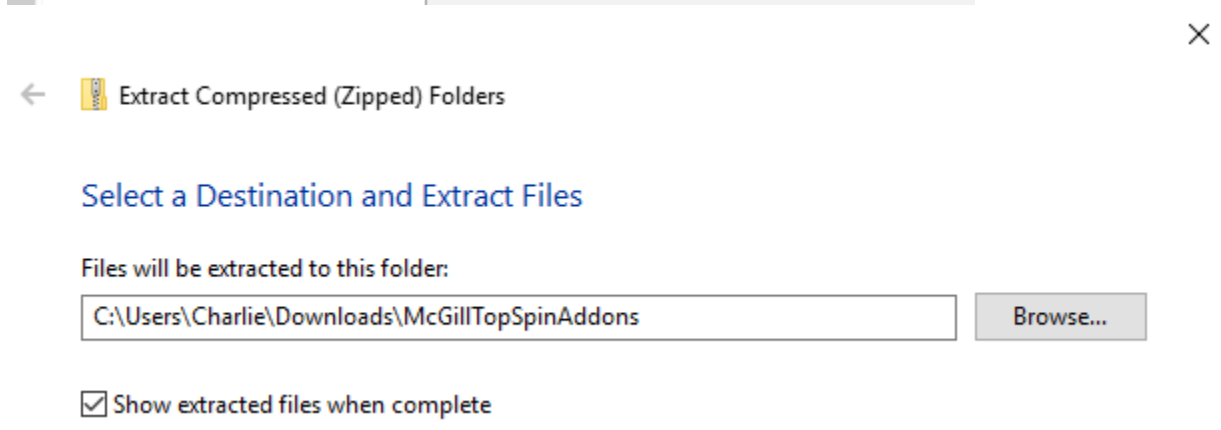


2. Open the folder containing McGillTopSpinAddons.zip and unzip it

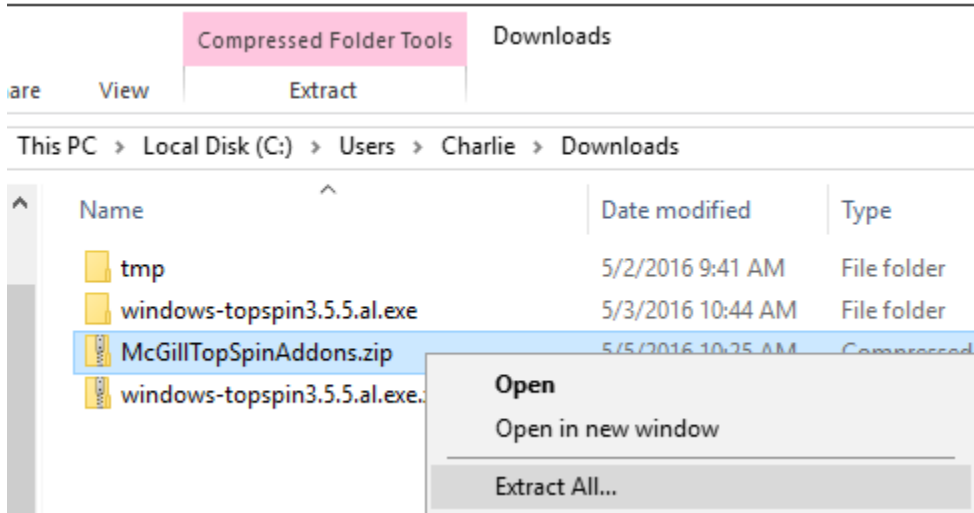


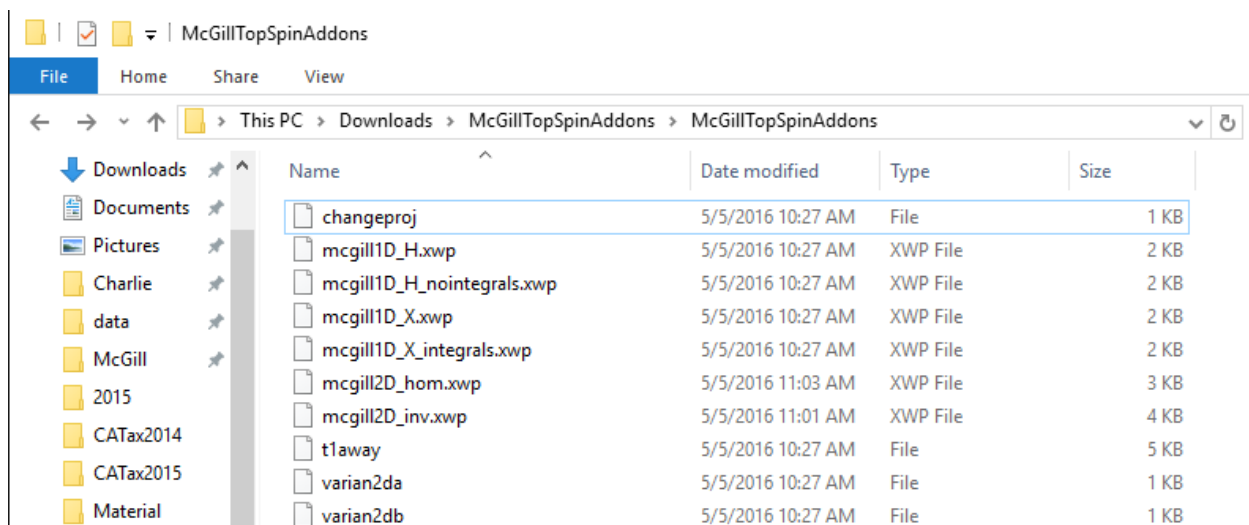


- Keep the Windows Explorer or Finder window showing the files open
- At the bottom of the TopSpin window, type `expl top`. This should open a Windows Explorer or Finder window showing the location of the TopSpin installation. If not, open Windows Explorer or the Finder and browse to the appropriate location (`c:\bruker\topspin4.0.5` or `/opt/topspin`)

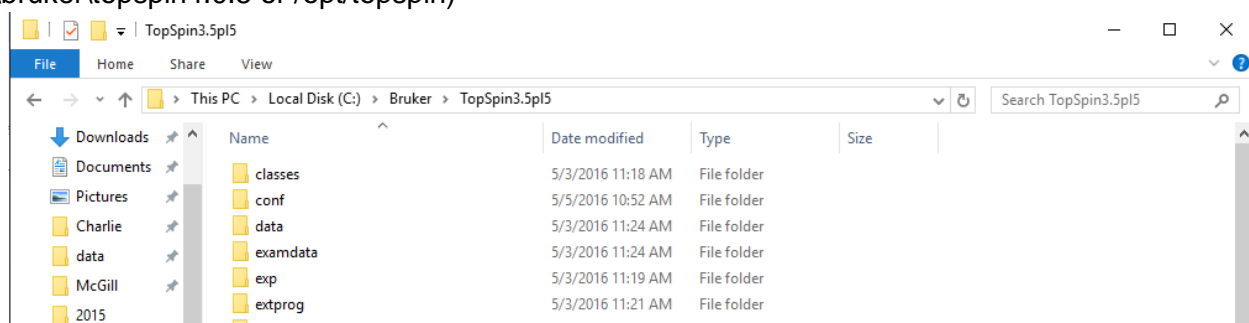


- From these, browse to the directory plot / layouts

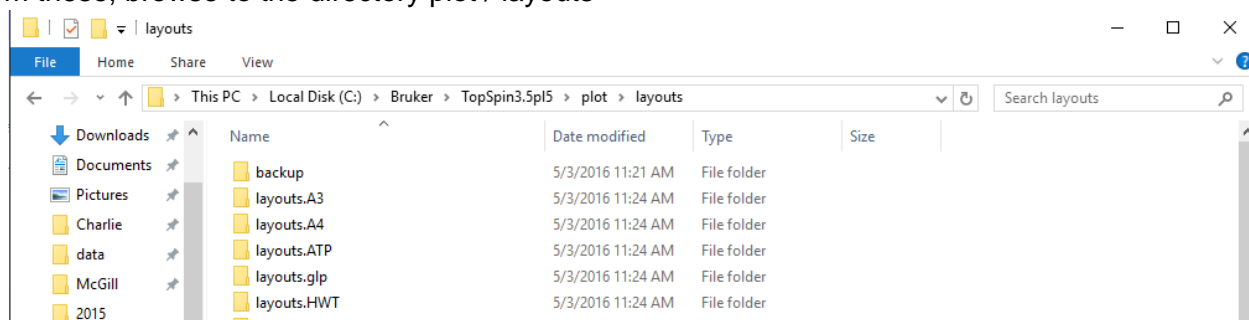

- Copy the files ending in `.xwp` from the zip file extracted in step 2 into this directory

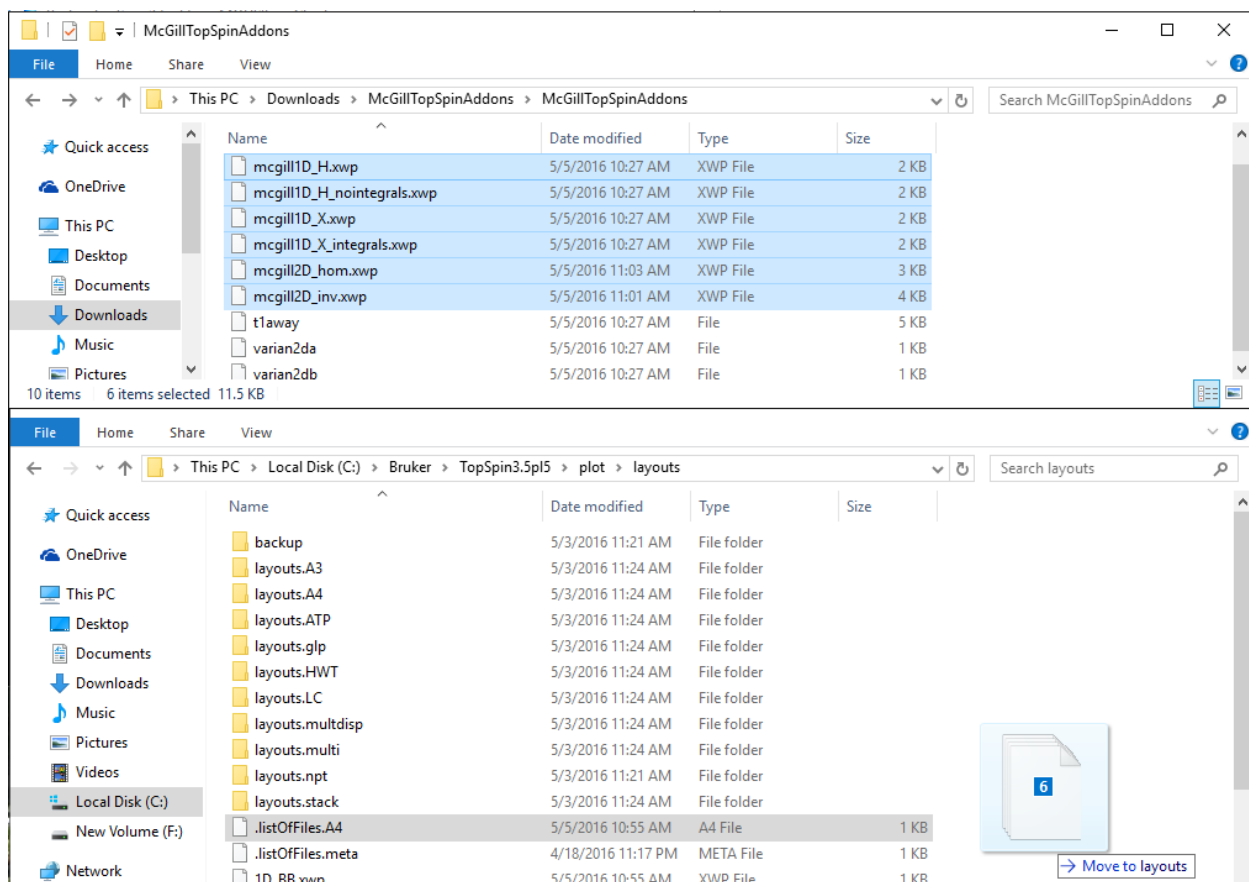

7. Next, go up to the TopSpin directory, and then browse to exp / stan / nmr / au / src / user

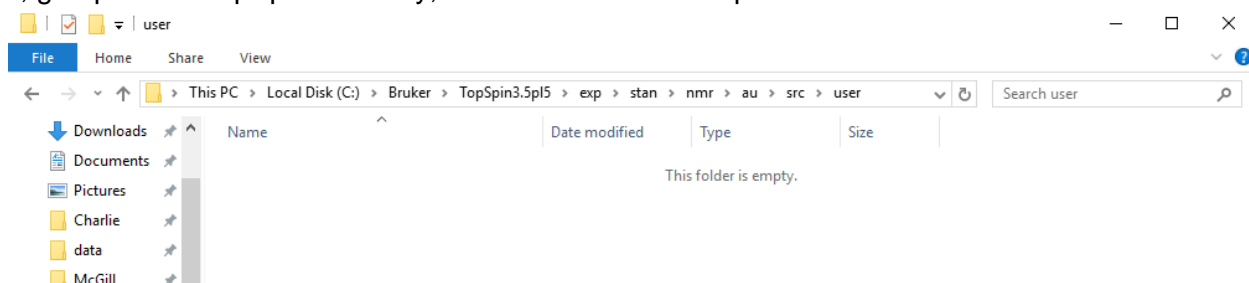

8. Copy all remaining files from the unzipped file into this directory

McGillTopSpinAddons

File Home Share View

This PC > Downloads > McGillTopSpinAddons > McGillTopSpinAddons


Name	Date modified	Type	Size
mcgill1D_H.xwp	5/5/2016 10:27 AM	XWP File	2 KB
mcgill1D_H_nointegrals.xwp	5/5/2016 10:27 AM	XWP File	2 KB
mcgill1D_X.xwp	5/5/2016 10:27 AM	XWP File	2 KB
mcgill1D_X_integrals.xwp	5/5/2016 10:27 AM	XWP File	2 KB
mcgill2D_hom.xwp	5/5/2016 11:03 AM	XWP File	3 KB
mcgill2D_inv.xwp	5/5/2016 11:01 AM	XWP File	4 KB
t1away	5/5/2016 10:27 AM	File	5 KB
varian2da	5/5/2016 10:27 AM	File	1 KB
varian2db	5/5/2016 10:27 AM	File	1 KB

10 items 4 items selected 5.84 KB

File Home Share View

This PC > Local Disk (C:) > Bruker > TopSpin3.5p15 > exp > stan > nmr > au > src > user

This folder is empty.

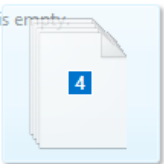

→ Move to user