

Users will now be fully autonomous for the use of the Nanodrop and Enspire. Please follow those instructions for your safety and the safety of others.

1. Please reserve your spot on this website (if you don't have a username and password, just contact me):

<https://mcgill.ca/c-bigneuro/core-facility/equipment>

2. Wear your PPE equipment (labcoat, gloves, mask)

3. Wash your hands before using the equipment or wear gloves that you disinfect with ethanol 70%.

4. Wash the equipment and bench with ethanol 70% (put some on a paper towel and wipe everything (even the keyboard and mouse)).

5. Turn on the equipment and log in into the computer (password: bluestar2020)

6. Carry on with your experiment.

7. Fill out the logbook (Nanodrop: Excell file on the computer, Enspire Plate Reader: Sheet on the reader - extra ones are in the sheet protector on the wall).

8. Turn off the equipment

9. Wash the equipment and bench with ethanol 70% (put some on a paper towel and wipe everything even the keyboard and mouse).

Thank you for your cooperation and be safe! If any questions or problems, please contact cbig at cbig.mni@mcgill.ca