

Instructions for Staining Polyacrylamide Gels

SDS-PAGE Mini Gels:

1. Wash gel 3 times for 5 min. each in 200 ml of ddH₂O per gel.
2. Remove all water from the staining container and add 50 ml Bio-Safe Coomassie Stain (or enough to completely cover gel). Gently shake for 1 hr. Protein bands will be visible within 20 min. and reach maximum intensity within 1 hr. Longer incubation will not increase background.
3. Rinse gel in 200 ml of ddH₂O for at least 30 min. Stained gels can be stored in water.

Note: Residual SDS in the gel may cause background staining and interfere with band intensity while the gel is in the stain. Rinsing the gel extensively in water after staining will remove background and allow proper visualization of the bands.

SDS-PAGE Gels:

For larger gels (16-20") increase volumes in above method 2 fold.

PEPTIDE Gels:

Fix gels in 40% methanol, 10% acetic acid for 30 min., follow with staining starting with Step 2 (above). Extend Step 3 (above) water wash to 2 hrs. Peptide bands will not be clearly visible until after the final water wash.

Ordering Information

Catalog #	Description
161-0786	Bio-Safe Coomassie Stain, 1L
161-0787	Bio-Safe Coomassie Stain, 5L

For technical service, contact your local Bio-Rad office, or in the US, call **1-8004BIORAD (1-800-424-6723)**

Bio-Rad Laboratories, Inc.
2000 Alfred Nobel Dr., Hercules, CA 94547
510-741-1000

4110120 Rev A